

Vion IMS QToF Shows Three Orders of Magnitude Linearity of Response for Metabolic Profiling

Adam King and Chris Hughes
 Waters Corporation, Milford, MA, USA



GOAL

To outline the quantification capabilities and linear response of the Vion IMS QToF for the analysis of human urine.

BACKGROUND

Metabolic profiling is a fast-emerging science which provides insight into changes in biochemical pathways in response to disease state or a toxicological insult. It provides insight not only into which metabolites have been up or down regulated, but the rate of these changes. Successful metabolomics methodologies involve the detection and quantification of hundreds or thousand of metabolites in biochemical fluids such as plasma and urine.

Accurate mass LC-MS has become a mainstay of metabolomic research due to its sensitivity, specificity, speed of analysis, and structural identification abilities. In order to provide reliable quantitative data, the mass spectrometry system must provide a linear response over the systemic concentration range of the metabolites.

The Vion™ IMS QToF is equipped with a novel two-stage ToF detector, QuanTof 2 Technology, allowing the detection of

The demonstration of Vion IMS QToF linearity of response for metabolic profiling in biomedical research.

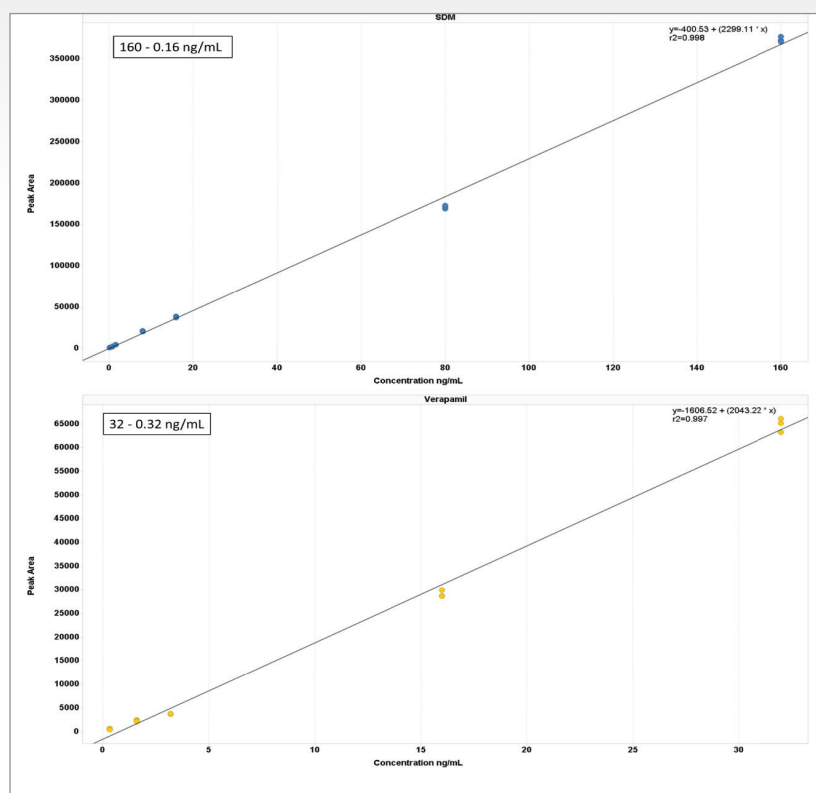


Figure 1. Calibration lines obtained for sulfadimethoxine (SDM) and verapamil in human urine.

metabolites over a wide dynamic range. Here we demonstrate the linear dynamic range of the Vion IMS QToF Mass Spectrometer for metabolomics analysis.

THE SOLUTION

The linear dynamic range of the Vion IMS QToF was evaluated in urine spiked with the Waters™ LC-MS test mix, containing seven non-endogenous compounds diluted to achieve a range over seven orders of magnitude. The urine samples were subject to protein precipitation with acetonitrile prior to analysis by UPLC-IMS-MS.

The samples were chromatographed on an ACQUITY™ UPLC™ BEH Amide, 1.7 μm, 2.1 x 100 mm Column (p/n: [186004849](#)), operated under HILIC gradient conditions over 10 min. The mobile phases consisted of 10 mM ammonium formate and formic acid in 95:5 water:acetonitrile (mobile phase A) and 5:95 water:acetonitrile (mobile phase B). The column effluent was monitored using a Vion IMS QTof operating in HDMS^E positive or negative mode with a collision energy ramp from 20–30 eV and a mass range of m/z 50–1200.

A typical calibration line obtained for the sulfadimethoxine and verapamil are shown below in Figure 1 and reserpine and sulfagaunidine in Figure 2. The data acquired shows linear response over four orders of magnitude for all the spiked compounds.

SUMMARY

Metabolic profiling provides a unique insight into biochemical changes as a result of disease or toxicity. The biological variation that can be seen when comparing diseased subjects to those of healthy controls produces a broad response from metabolites often to the extremes of instrument capabilities. The UPLC-MS-MS system showed a linear dynamic range over three orders of magnitude for small molecules in urine, making the system ideal for the semi-quantitative measurement and identification of metabolites in biological fluids.

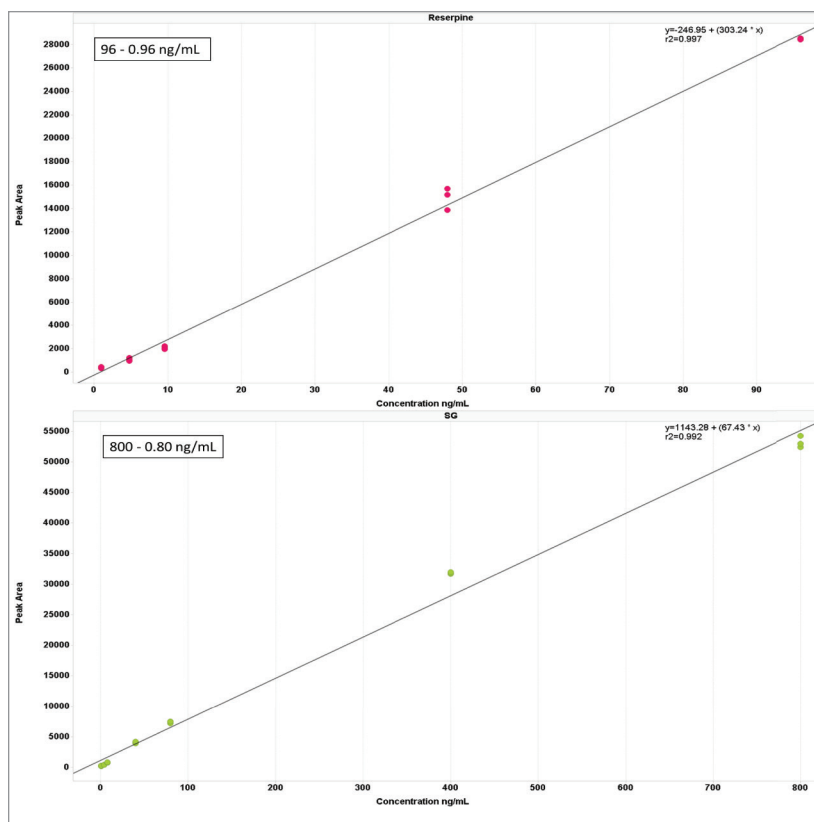


Figure 2. Calibration line obtained for reserpine and sulfagaunidine (SG) in human urine.

For research use only. Not for use in diagnostic procedures.

Waters

THE SCIENCE OF WHAT'S POSSIBLE.™

Waters, The Science of What's Possible, ACQUITY, UPLC, and Vion are trademarks of Waters Corporation. All other trademarks are the property of their respective owners.

©2019 Waters Corporation. Produced in the U.S.A. September 2019 720006643EN TC-PDF

Waters Corporation
34 Maple Street
Milford, MA 01757 U.S.A.
T: 1 508 478 2000
F: 1 508 872 1990
www.waters.com