

MassHunter VistaFlux Ver 1.0 Site Preparation Checklist

Thank you for purchasing an Agilent instrument. To get you started and to assure a successful and timely installation of your Agilent **software**, please refer to this specification or set of requirements.

Correct site preparation is the key first step in ensuring that your instruments and software systems operate reliably over an extended lifetime. This document is an **information guide and checklist** prepared for you that outlines the supplies, consumables, space and utility requirements for your equipment.

Customer Responsibilities

Make sure your site meets the following prior specifications before the installation date. For details, see specific sections within this checklist, including:

- The computing environment and the necessary space is made available
- The number & location of electrical outlets for your computer systems and peripherals are planned
- That your site meets the software, hardware and networking specifications below
- Locate your sales order information, software authorization codes and/or software licenses/certificates
- The necessary software media, disks etc are available including upgrade/update disks
- That a suitable backup solution is identified for your software
- Availability of a system/network administrator as needed to connect to your intranet
- Please consult Other Requirements section below for other product-specific information.


If Agilent is delivering installation and familiarization services, users of the instrument should be present throughout these services; otherwise, they will miss important operational, maintenance and safety information.

- Connection to the internet is necessary to complete the installation of this software product.

Important Customer Information

1. If you have questions or problems in providing anything described as a Customer Responsibility above, please contact your local Agilent or partner support/service organization for assistance prior to delivery. In addition, Agilent and/or its partners reserve the right to reschedule the installation dependent upon the readiness of your laboratory.
2. Should your site not be ready for whatever reasons, please contact Agilent as soon as possible to re-arrange any services that have been purchased.
3. Other optional services such as additional training, and consultation for user-specific applications may also be provided at the time of installation when ordered with the system, but should be contracted separately.

**MassHunter VistaFlux Ver 1.0
Site Preparation Checklist**

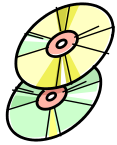

**Software Specifications for
Workstations, Clients, and Servers**

Special Notes

1. Author to add special considerations or notes.

Specification Description	Minimum	Recommended (if applicable)	Comments (if applicable)
Operating system type(s), versions	Microsoft Windows 7 Professional SP1 64-Bit		
O/S .NET & other add-ons	Microsoft® NET 3.5 Service Pack 1		
Language settings/compatibility	English		
Regional settings/compatibility	English		
Account settings/privileges	Administrator privileges required for installation.		
Specific drivers	Not Applicable		



**Computer Hardware Specifications for
Workstations, Clients, and Servers**

Special Notes

1. Author to add special considerations or notes.

Specification Description	Minimum	Recommended (if applicable)	Comments (if applicable)
Processor type & speed	Intel Core i5 > 2.0 GHz	4 Core >3.0 GHz	
Memory	16 Gigabyte	24- 32 Gigabyte	24 to 32 GB RAM is recommended for improved data processing in Profinder B.08 or higher.
Internal Storage/devices/media	1 Terabyte		Depends on size of projects and datasets
External Storage/devices/media	Not Applicable		
Video devices	SVGA or better graphics adapter and monitor		

**MassHunter VistaFlux Ver 1.0
Site Preparation Checklist**

	(minimum resolution 1280 x 1024), True Colors		
Audio devices	Not Applicable		
Printing devices	Not Applicable		
Pointing devices	Not Applicable		


Networking Specifications
Special Notes

- Check as not applicable.

Specification Description	Minimum	Recommended (if applicable)	Comments (if applicable)
Network type, bandwidth, speed, protocol etc	Not Applicable		
Additional network or instrument devices/cards requirements	Not Applicable		


Other Requirements

Author to add other guidance, notes and/or diagrams that are needed by the customer which have not been mentioned above such as:

- Important web links
Microsoft Hardware Compatibility Lists
<http://www.microsoft.com/whdc/hcl/default.msp>

Links to specific O/S fixes, updates needed
<http://support.microsoft.com/>
- Additional or supporting applications/tools, diagnostics or add-ons needed or recommended.
- Additional set up guides or manuals to be referenced.
- Additional info regarding required firmware for instruments.
- Additional upgrade and/or migration preparation instructions.

**MassHunter VistaFlux Ver 1.0
Site Preparation Checklist**

Important Customer Web Links

- For additional information about our solutions, please visit our web site at <http://www.agilent.com/en-us/solutions>
- Need to get information on your product?
Literature Library - <http://www.agilent.com/chem/library>
- Need to know more?
Customer Education - <http://www.agilent.com/chem/education>
- Need technical support, FAQs? - <http://www.agilent.com/chem/techsupp>
- Need supplies? - <http://www.agilent.com/chem/supplies>

Document part number: G4492-90005