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## Dear Customer,

Thank you for purchasing a product from Agilent Technologies. Before you start to use your product, please read the corresponding documentation provided on this USB flash drive. It provides an extensive collection of user manuals, safety information and technical reference for individual instruments.

## General Safety

The following general safety precautions must be observed during all phases of operation of this instrument. Failure to comply with these precautions or with specific warnings or operating instructions in the user manuals violates safety standards of design, manufacture, and intended use of the instrument. Agilent Technologies assumes no liability for the customer's failure to comply with these requirements. User manuals are provided with your instrument on this USB flash drive and/or in printed form. Manuals may also be available on the Web.

Go to **www.agilent.com** and type in your product number in the Search field at the top of the page.

## General

Do not use this product in any manner not specified by the manufacturer. The protective features of this product may be impaired if it is used in a manner not specified in the operation instructions.

## Before Applying Power

Verify that all safety precautions are taken according to the user manual of the individual instrument. Verify that the voltage range and frequency of your power distribution matches to the power specification of the individual instrument. Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations. Make all connections to the unit before applying power. Note the instrument's external markings described under "Safety Symbols" in the Appendix of the corresponding user manual.

## Ground the Instrument

If your product is provided with a grounding type power plug, the instrument chassis and cover must be connected to an electrical ground to minimize shock hazard. The ground pin must be firmly connected to an electrical ground (safety ground) terminal at the power outlet. Any interruption of the protective (grounding) conductor or disconnection of the protective earth terminal will cause a potential shock hazard that could result in personal injury.

## Instrument Connection

Any other instruments connected to this instrument shall be approved to a suitable safety standard and must include reinforced insulation from the mains.

## Do Not Operate in an Explosive Atmosphere

Do not operate the instrument in the presence of flammable gases or fumes.

## Do Not Remove the Instrument Cover

Only Agilent authorized personnel are allowed to remove instrument covers.

## Do Not Modify the Instrument

Do not install substitute parts or perform any unauthorized modification to the product. Return the product to Agilent or consult an Agilent Sales and Service Office for service and repair to ensure that safety features are maintained.

## In Case of Damage

Instruments that appear damaged or defective should be made inoperative and secured against unintended operation until they can be repaired by qualified service personnel.

## Toxic and hazardous solvents and flammable liquids

The handling of solvents and reagents can hold health risks. When working with solvents observe appropriate safety procedures (for example, goggles, safety gloves and protective clothing) as described in the material safety data sheet supplied by the solvent vendor, especially when toxic or hazardous solvents and flammable liquids are used.

Do not use solvents with an auto-ignition temperature below 200°C (392°F).

Do not use solvents with a boiling point below 56°C (133°F).

Avoid high vapor concentrations.

Always keep the solvent temperature at least 25 K below the boiling point of the solvent used.

Do not use solvents of ignition Class IIC according IEC 60079-20-1 (for example, carbon disulfide).

Reduce the volume of substances to the minimum required.

Ground the waste container.

Regularly check the filling level of the waste container. The residual free volume in the waste container must be large enough to collect the waste liquid.

To achieve maximal safety, regularly check the tubing for correct installation.

Note that for some Agilent products, precautions and limits apply that may differ from the specifications above. Always follow the relevant instructions shipped with a specific product.

## Safety Notices

### CAUTION

A **CAUTION** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in damage to the product or loss of important data. Do not proceed beyond a **CAUTION** notice until the indicated conditions are fully understood and met.

### WARNING

A **WARNING** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in personal injury or death. Do not proceed beyond a **WARNING** notice until the indicated conditions are fully understood and met.

## Safety Symbols

Symbol	Description
 Caution	The apparatus is marked with this symbol when the user should refer to the instruction manual in order to protect risk of harm to the operator and to protect the apparatus against damage.
 electrical hazard	Indicates dangerous voltages.
 pinching or crushing	Indicates a pinching or crushing hazard.
 piercing or cutting	Indicates a piercing or cutting hazard.
 Magnetic field	Magnets produce a far-reaching, strong magnetic field. They could damage TVs and laptops, computer harddrives, credit and ATM cards, data storage media, mechanical watches, hearing aids and speakers. Keep magnets at least 25 mm away from devices and objects that could be damaged by strong magnetic fields.
 pacemaker	Magnets could affect the functioning of pacemakers and implanted heart defibrillators. A pacemaker could switch into test mode and cause illness. A heart defibrillator may stop working. If you wear these devices keep at least 55 mm distance to magnets. Warn others who wear these devices from getting too close to magnets.
 hot surface	The apparatus is marked with this symbol when hot surfaces are available and the user should not touch it when heated up.
 flammable	For Sampler Cooler which uses flammable refrigerant consult Agilent Information Center / User Manual before attempting to install or service this equipment. All safety precautions must be followed.
	Confirms that a manufactured product complies with all applicable European Community directives. The European Declaration of Conformity is available at: <a href="http://regulations.corporate.agilent.com/DoC/search.htm">http://regulations.corporate.agilent.com/DoC/search.htm</a>
	Manufacturing date.
	Power symbol indicates ON/Off. The apparatus is not completely disconnected from the mains supply when the power switch is in the Off position.
	Cooling unit is designed as vapor-compression refrigeration system. Contains fluorinated greenhouse gas (refrigerant) according to the Kyoto protocol.
	Indicates a protected ground terminal.
 Laser	Internal Laser beam label located on light source.

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© Agilent Technologies, Inc. 2014–2023