


ISOMIST™ – BLUETOOTH SET UP GUIDE

Windows® XP / Windows® 7 / Windows® 8

Before you start: The IsoMist application should already be installed on your PC. Ensure the USB cable is not connected to the IsoMist.



1

Plug the Bluetooth dongle into a free USB port on your PC. Your computer will automatically detect and install the drivers for the Bluetooth dongle.

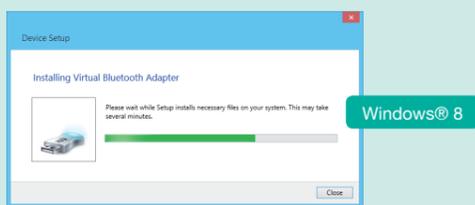

2




Wait until the Bluetooth dongle has been installed.

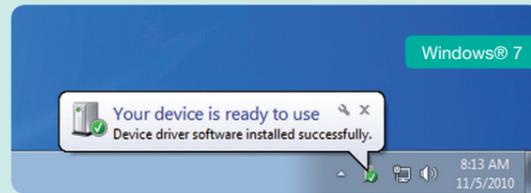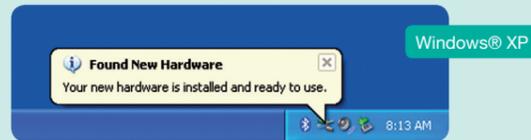

Windows® 8, the drivers have been installed when the windows close.




3

Plug in and turn on the IsoMist, ensuring that it is not connected through the USB cable. Within one minute a Bluetooth balloon pop-up will appear. Click this balloon pop-up.

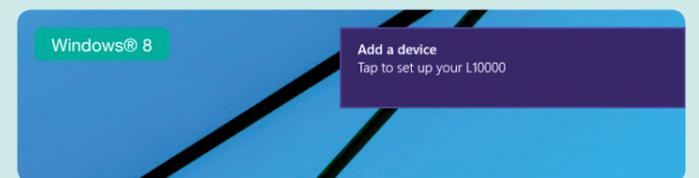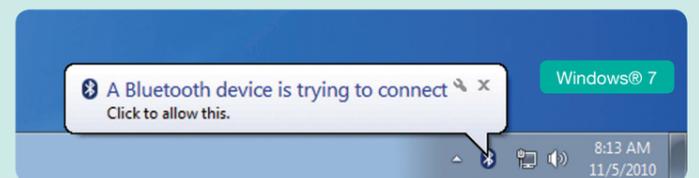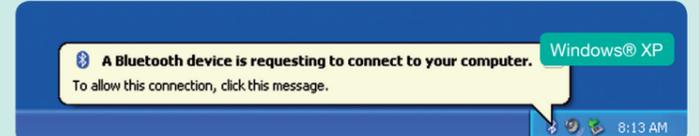



4

Enter the passkey / pairing code for the IsoMist in the **Add Device** window that appears, and click **Next**. The passkey is the same as the serial number of your IsoMist. This can be found on the **Add Device** window itself, next to a graphical icon. It can also be found on the underside of your IsoMist.

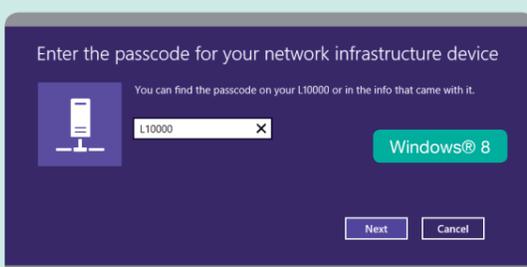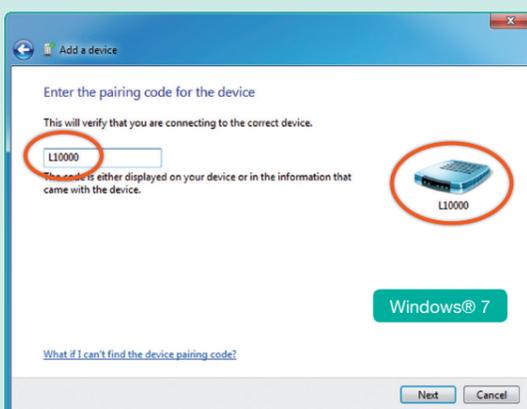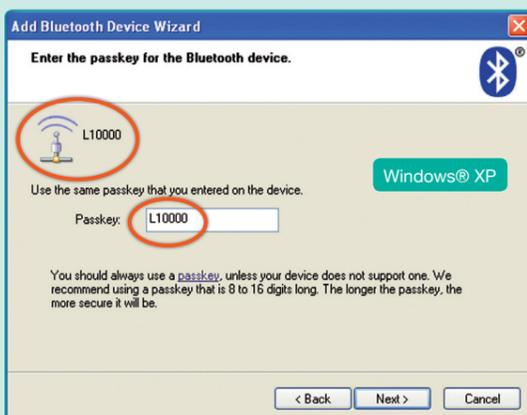

5

The following window confirms that the IsoMist has been paired to your PC. Click **Close / Finish**.

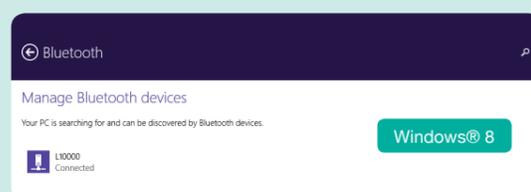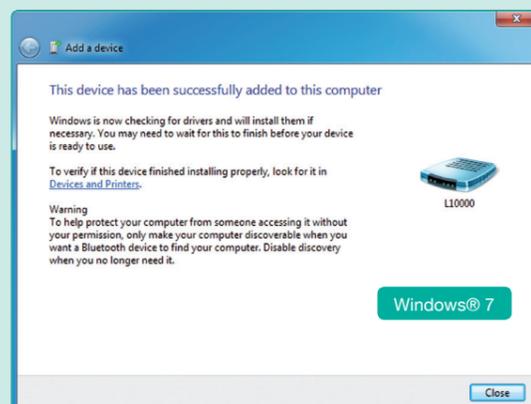


6

Launch the IsoMist software from your desktop, or the Start menu. Once the IsoMist software is open, click the **Power** button to connect to the IsoMist.



7

The **Connect** button in the IsoMist software will now change to a **Bluetooth** button, to verify that the IsoMist is connected and operational via Bluetooth communications.

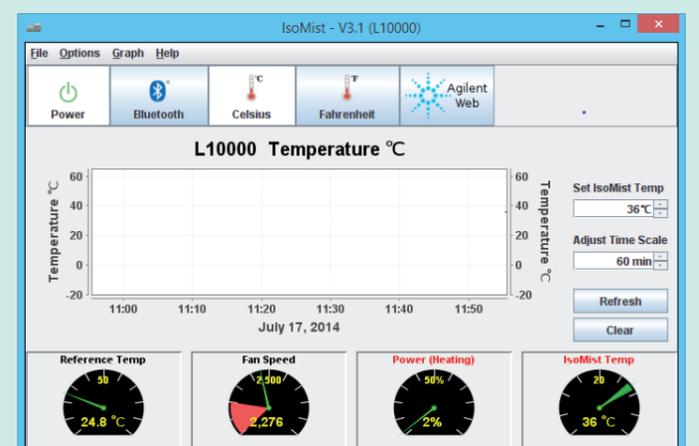


ISOMIST™ – SOFTWARE INSTALLATION GUIDE

Windows® XP / Windows® 7 / Windows® 8


Before you start: Ensure your user account has Administrative privileges

START


1. Exit all other programs you may have running on your PC.
2. Insert the IsoMist software installation memory stick (supplied with IsoMist pack) into a free USB port on your PC.
3. Locate and run “Setup.exe” on the memory stick.

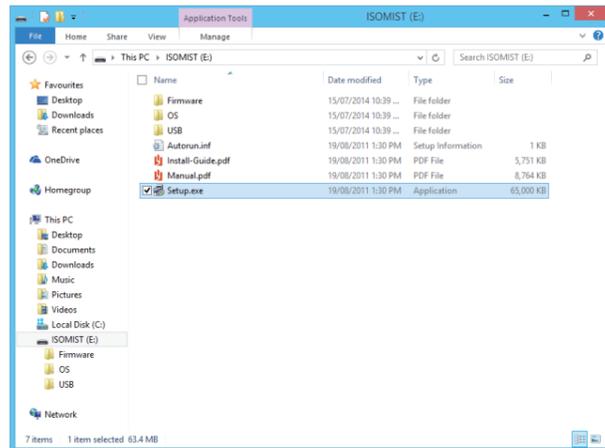


4. If the following dialogue appears, click **Yes**.

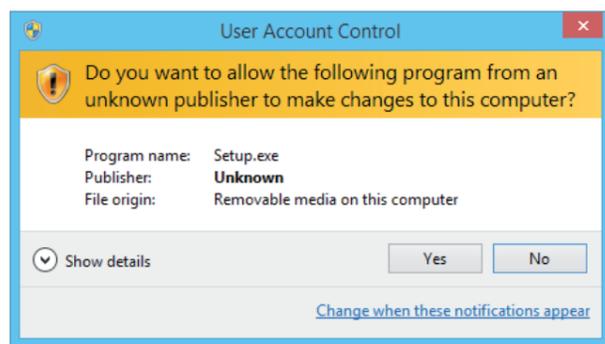


5. Click **Yes**.

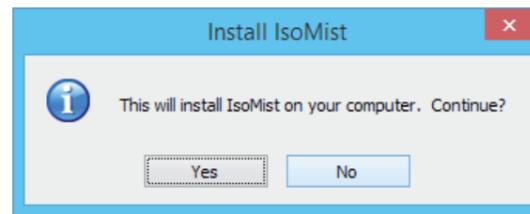


6. Click **Next**.

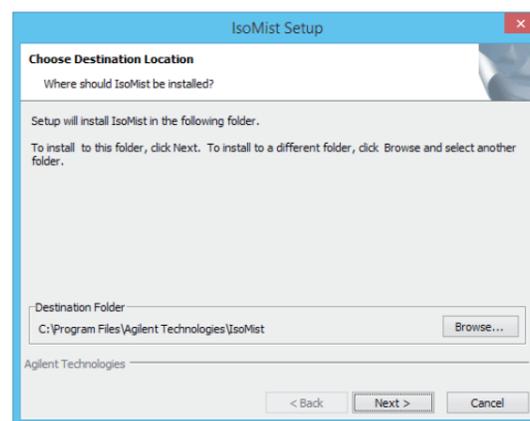


7. Click **Next**.

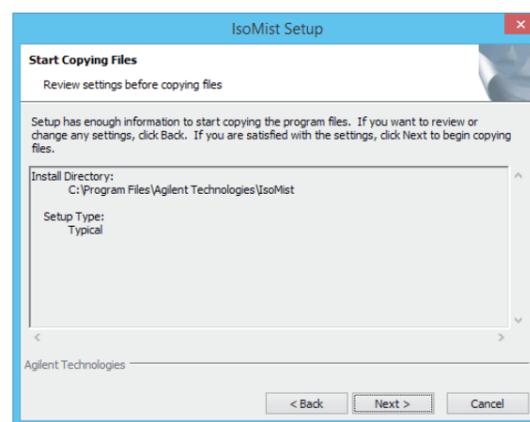

8. When the device driver wizard pops up, click **Next**

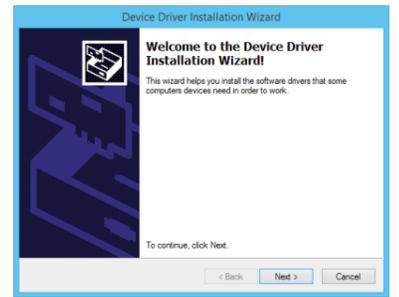

9. Windows® 7: Make sure **Always trust software from “Belcarra Technologies (2005) Corp.”** is checked, and click **Install**.

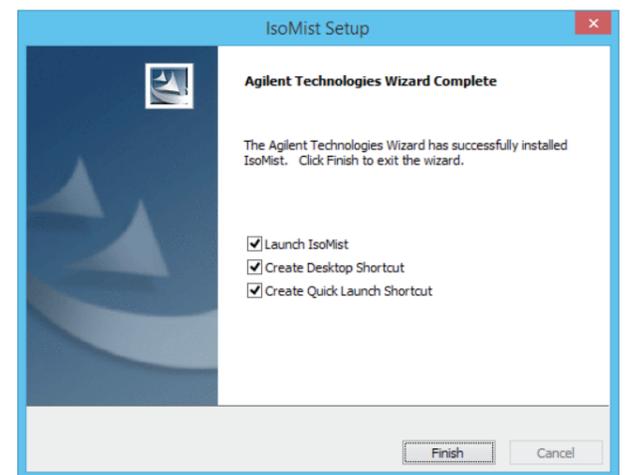

10. The first time IsoMist is plugged into a USB port, it will need to activate the USB drivers.



This is done automatically in Windows® 7, and the following message will appear to indicate successful activation.




For Windows® XP activation, please see the following steps:

After IsoMist is plugged into a USB port for the first time, the USB driver must first be activated.

1. When the following window appears, select **No, not this time**, and click **Next**.
2. When the following window appears, select **Install the software automatically**, and click **Next**.
3. If the following window appears, click **Continue Anyway**.
4. When the following window appears, click **Finish**.

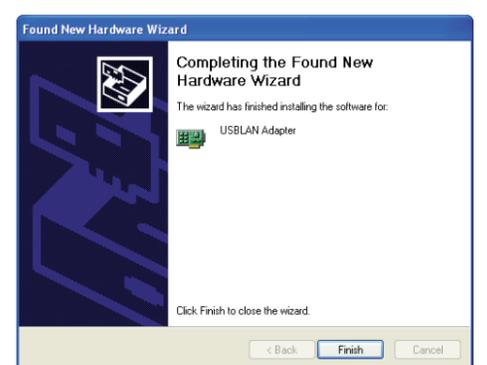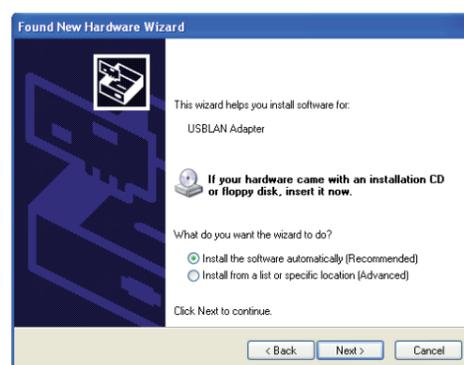

5. Windows® XP driver activation is now complete. The following message will appear to indicate successful activation.

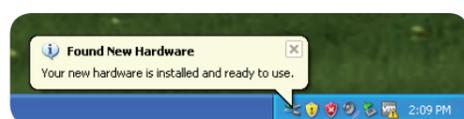

Agilent Technologies