

Generic Method for the Extraction of Quaternary Amines and Polybasic Drugs from Biological Fluids Using ISOLUTE® HCX-Q Mixed-Mode SPE Columns and 96-well Plates

Introduction

The use of ISOLUTE® HCX mixed-mode sorbents is widely accepted for providing high purity extracts of basic drugs from biological fluids. The dual retention mechanism, provided by non-polar and strong cation exchange functional groups, allows for a rigorous interference elution step resulting in low levels of co-extracted materials, thus minimizing the risk of ion suppression effects in LC-MS applications. This approach is suitable for drugs that can be eluted under basic conditions. However, quaternary amines and polybasic drugs cannot be eluted under these conditions. ISOLUTE® HCX-Q, containing non-polar and weak cation exchange functional groups, was designed to allow elution of these compounds under acidic elution conditions.

The ISOLUTE HCX-Q sorbent utilizes a combination of weak cation exchange and C8 non-polar retention mechanisms. The analyte is initially retained by a non-polar (hydrophobic) retention mechanism, which is unaffected by high or variable ionic strength of the sample matrix. The pH control and salt removal at the sample application stage ensures that the polybasic drug or quaternary amine is retained by a robust cation exchange retention mechanism. The two retention mechanisms are shown in Figure 1. An interference elution step containing both aqueous buffer and a water miscible organic solvent (e.g. methanol) is used to remove many non-polar interferences. The quaternary amines are eluted by breaking both non-polar and ionic interactions, providing recovery of the analyte.

Extraction Procedure

This procedure utilizes ISOLUTE® HCX-Q 25 mg/1 mL column or 96-well SPE plate configurations, optimized for the elution of basic drugs in very low volumes. Solvent volumes should be scaled up or down as necessary for other column configurations.

Sample Pre-treatment

Dilute the 100 µL biological fluid sample with 0.05 M ammonium acetate buffer, pH 7 (1:4, v/v). Mix thoroughly.

Column Conditioning and Equilibration

Condition the column with methanol (1 mL) followed by 0.05 M ammonium acetate buffer, pH 7 (500 µL).

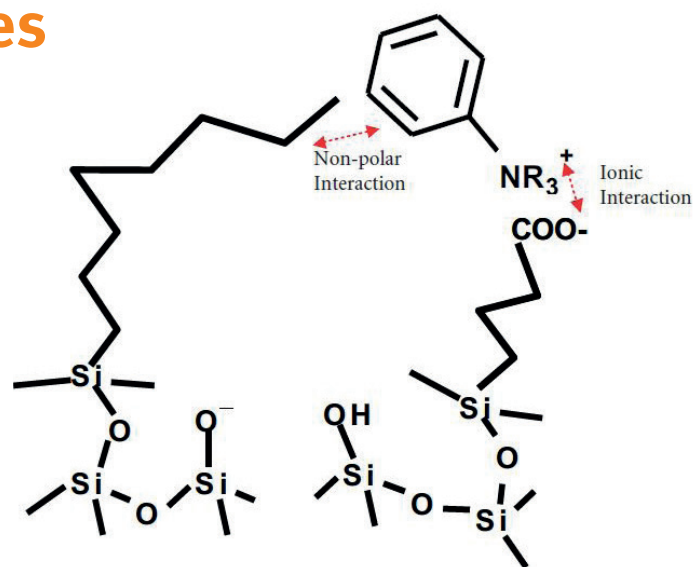


Figure 1. Multiple interactions on ISOLUTE® HCX-Q.

Sample Application

Apply the sample (500 µL) at a flow rate of 1–2 mL/minute.

Interference Elution

Rinse the column with methanol/0.05 M ammonium acetate buffer, pH 7.0 (20:80, v/v, 500 µL). The buffer ensures that the sorbent surface remains negatively charged during the interference elution step.

Increase and optimize the proportion of methanol to improve extract cleanliness, checking for analyte breakthrough.

Analyte Elution

Elute analytes with 0.2 M monochloroacetic acid in methanol (2 x 125 µL).

Analyte elution is achieved by simultaneously suppressing ionization of the sorbent carboxylic acid functional groups and eliminating non-polar interactions through the use of the acidified methanol elution solvent. The use of higher concentration acid in the elution solvent may be necessary for some analytes.

If the final analysis technique is GC, evaporate the elution solvent to dryness and derivatize the analyte(s) using a suitable derivatization agent.

This Technical Note is based on extraction conditions developed by Lindegardh et al¹ using a suite of antimalarial quaternary amine drugs.

References

¹Automated mixed-mode solid-phase extraction for simultaneous determination of atovaquone, proguanil and metabolites in plasma by HPLC.

Lindegardh, N; Bergqvist, Y. Department of Analytical Chemistry, Uppsala University, Sweden.

Presented at Analytical Days 2002, Sweden.

Ordering Information

Part Number	Description	Quantity
ISOLUTE® HCX-Q SPE Columns		
986-0002-A	ISOLUTE® HCX-Q 25 mg/1 mL columns	100
986-0010-A	ISOLUTE® HCX-Q 100 mg/1 mL columns	100
ISOLUTE-96 fixed well 96-well plates for high throughput		
986-0025-P01	ISOLUTE®-96 HCX-Q 25 mg Fixed well plate	1
986-0100-P01	ISOLUTE®-96 HCX-Q 100 mg Fixed well plate	1

EUROPE

Main Office: +46 18 565900
Toll Free: +800 18 565710
Fax: +46 18 591922
Order Tel: +46 18 565710
Order Fax: +46 18 565705
order@biotage.com
Support Tel: +46 18 56 59 11
Support Fax: +46 18 56 57 11
eu-1-pointsupport@biotage.com

NORTH & LATIN AMERICA

Main Office: +1 704 654 4900
Toll Free: +1 800 446 4752
Fax: +1 704 654 4917
Order Tel: +1 704 654 4900
Order Fax: +1 434 296 8217
ordermailbox@biotage.com
Support Tel: +1 800 446 4752
Outside US: +1 704 654 4900
us-1-pointsupport@biotage.com

JAPAN

Tel: +81 3 5627 3123
Fax: +81 3 5627 3121
jp_order@biotage.com
jp-1-pointsupport@biotage.com

CHINA

Tel: +86 21 68162810
Fax: +86 21 68162829
cn_order@biotage.com
cn-1-pointsupport@biotage.com

KOREA

Tel: +82 31 706 8500
Fax: +82 31 706 8510
korea_info@biotage.com
kr-1-pointsupport@biotage.com

INDIA

Tel: +91 11 45653772
india@biotage.com

Distributors in other regions
are listed on www.biotage.com

Literature Number: TN129.V.1

© 2020 Biotage. All rights reserved. No material may be reproduced or published without the written permission of Biotage. Information in this document is subject to change without notice and does not represent any commitment from Biotage. E&OE. A list of all trademarks owned by Biotage AB is available at www.biotage.com/legal. Other product and company names mentioned herein may be trademarks or registered trademarks and/or service marks of their respective owners, and are used only for explanation and to the owners' benefit, without intent to infringe.