

## Residual Solvents Analysis in Pharmaceuticals by HS-GC-FID with Newly Added Compounds – USP <467> Procedure A

Elgin Ting<sup>1</sup>, Cynthia Lahey<sup>1</sup>  
1 Shimadzu (Asia Pacific) Pte Ltd.

### User Benefits

- ◆ Good separation was achieved with newly added compounds
- ◆ Cost effective when using nitrogen as carrier gas instead of helium

### Introduction

United States Pharmacopeia (USP) <467> Residual Solvents is a compendium that defines acceptable amounts of residual solvents in pharmaceuticals and provides procedures to analyze residual solvents [1]. The recommended limit or classification of the solvent in USP <467> would be updated when new safety data on more solvents become available [2]. Methyl isobutyl ketone was reclassified as a Class 2 solvent in USP <467> in 2020 (previously a Class 3 solvent). Meanwhile, tertiary butyl alcohol and cyclopentyl methyl ether were proposed to be classified as Class 2 residual solvents in 2021 [1][2]. Here, we describe headspace-gas chromatography (HS-GC) method using nitrogen carrier gas to analyse Class 1 and Class 2 solvents, following USP<467> Procedure A, updated with 3 new solvents, namely cyclopentyl methyl ether, tertiary butyl alcohol and methyl isobutyl ketone.

### Measurement Conditions and Samples

HS-20 NX headspace autosampler paired with Nexis GC-2030 (Shimadzu Corporation, Japan) were used in this work. The analytical conditions following Procedure A under water-soluble article section in USP<467> are shown in Table 1.

Certified USP<467> Class 1 and 2 mixed standard solutions (without cyclopentyl methyl ether, tertiary butyl alcohol and methyl isobutyl ketone) were purchased from Restek Corporation, USA. The standards were prepared according to USP<467> Procedure A. Cyclopentyl methyl ether, tertiary butyl alcohol and methyl isobutyl ketone were purchased from Tokyo Chemical Industry (TCI). A mixture containing these 3 compounds was prepared in dimethyl sulfoxide at these concentrations: 7,500 µg/ml (cyclopentyl methyl ether), 17,500 µg/ml (tertiary butyl alcohol) and 22,500 µg/ml (methyl isobutyl ketone). Then, this mixture and Class 2A standard solution were mixed and further diluted following USP <467> Procedure.

### Results

#### Class 1 Standard


Class 1 result is displayed in Table 2. The average S/N value (n=7) for 1,1,1-trichloroethane (peak 2) is 114, which is much greater than the requirement stated in USP<467> (i.e., S/N ratio is not less than 5). Carbon tetrachloride (peak 3), which sensitivity is the lowest among the Class 1 solvents, has an average S/N value of 8. The repeatability of peak areas, %RSD (n=7) obtained from the 5 solvents ranges from 3.6% to 5.3%. These results indicate that this HS-GC method using nitrogen carrier gas can achieve the sensitivity requirement as stated in the USP<467> Procedure A for Class 1 Standards.

**Table 1:** HS-GC analytical conditions for residual solvent analysis following USP <467>

Instruments and Column information	
GC-FID	Nexis GC-2030
Auto Injector	HS-20 NX
Column	SH-I-624SIL MS 30 m x 0.32 mm ID x 1.80 µm df
Detector	FID-2030 Flame Ionization Detector
HS parameter	
Oven Temperature	80 °C
Sample Line Temperature	110 °C
Transfer Line Temperature	120 °C
Injection Time	1 min
Pressurizing Gas Pressure	75 kPa
Equilibrating Time	60 min
Shaking Level	2
GC-FID parameter	
Injection Mode	Split mode Split ratio 5
Carrier Gas	Nitrogen
Gas Flow Condition	Constant linear velocity mode Linear velocity 35 cm/s
Oven Temperature Programming	40 °C (20 min) → 10 °C/min to 240 °C (20 min)
Detector Temperature	250 °C
Hydrogen Flow	32 mL/min
Synthetic Air Flow	200 mL/min
Make-up Gas Flow	24 mL/min

**Table 2:** Peak area repeatability and average signal to noise ratio (S/N) for Class 1 Standard

Peak No.	Solvent	%RSD of peak area (n=7)	Average S/N ratio (n=7)
1	1,1-Dichloroethene	4.6	118
2	1,1,1-Trichloroethane	3.6	114
3	Carbon tetrachloride	5.3	8
4	Benzene	4.0	146
5	1,2-Dichloroethane	4.1	56

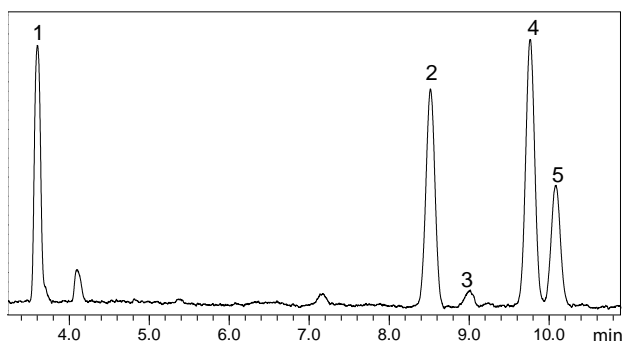

**Figure 1:** HS-GC-FID chromatogram of Class 1 Standard following Procedure A in USP<467>. Peak labelling refers to Table 2.

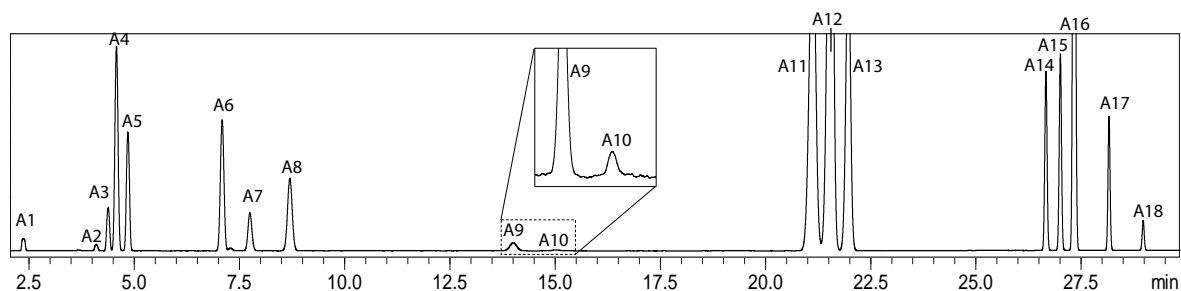


Figure 2: Chromatogram of Class 2A Standard analysed using Procedure A. Refer to Table 3 for the peak labeling.

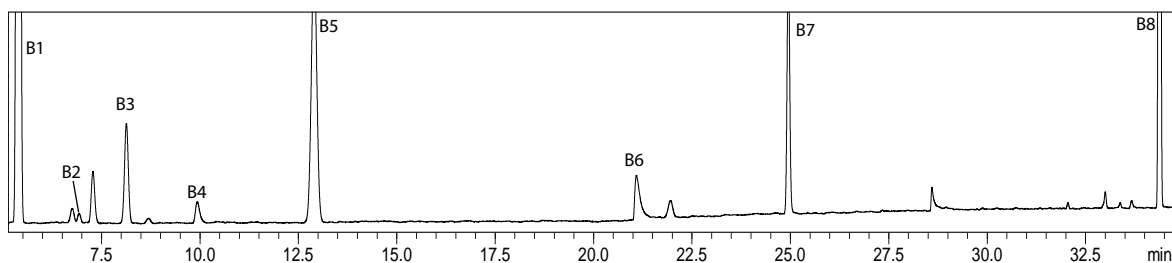


Figure 3: Chromatogram of Class 2B Standard analysed using Procedure A. Refer to Table 3 for the peak labeling.

### Class 2 Standard

Class 2 Standard consists of two groups, Class 2A and Class 2B. The chromatograms are shown in Figure 2 and 3, respectively. The repeatability results of Class 2A and Class 2B are compiled into Table 3. The peak area %RSD (n=7) values obtained for all the solvents ranges from 0.6% to 6.7%.

Acetonitrile and methylene chloride were well separated as shown in Figure 2. The specific resolution ( $R_s$ ) between acetonitrile and methylene chloride obtained was 2.2. This meets the stated criteria of USP<467> that  $R_s$  of these 2 compounds must not be less than 1. Even with three new compounds added into Class 2A standard mixture, all the peaks were well separated (Figure 2).

### Conclusion

HS-20 NX with Nexis GC-2030 with nitrogen carrier gas successfully achieved the required sensitivity (S/N) and peak resolution stated in the criteria of USP<467> Procedure A. Good separation of Class 2A solvents was maintained even with the addition of three new compounds, namely cyclopentyl methyl ether, tertiary butyl alcohol and methyl isobutyl ketone.

### References

1. The United States Pharmacopeia, USP <467> RESIDUAL SOLVENTS.
2. The International Council for Harmonisation of Technical Requirements for Pharmaceuticals for Human Use 2021, IMPURITIES: GUIDELINE FOR RESIDUAL SOLVENTS Q3C(R8), accessed 15 December 2021 <[https://database.ich.org/sites/default/files/ICH\\_Q3C-R8\\_Guideline\\_Step4\\_2021\\_0422\\_1.pdf](https://database.ich.org/sites/default/files/ICH_Q3C-R8_Guideline_Step4_2021_0422_1.pdf)>

Table 3: Peak area repeatability (n=7) for Class 2A and 2B Standards

Class 2A standard		
No.	Compound	%RSD (n=7) of peak area
A1	Methanol	0.6
A2	Acetonitrile	0.7
A3	Methylene chloride	3.2
A4	Tertiary Butyl Alcohol	0.6
A5	trans-1,2-Dichloroethylene	3.8
A6	Cis-1,2-Dichloroethylene	3.3
A7	Tetrahydrofuran	0.7
A8	Cyclohexane	3.8
A9	Methylcyclohexane	4.0
A10	1,4-Dioxane	6.7
A11	Cyclopentyl Methyl Ether	1.7
A12	Methylisobutyl Ketone	1.2
A13	Toluene	3.2
A14	Chlorobenzene	3.1
A15	EthylBenzene	3.2
A16	m,p-Xylene	3.1
A17	o-xylene	3.0
A18	Cumene	3.2
Class 2B standard		
No.	Compound	%RSD (n=7) of peak area
B1	Hexane	6.7
B2	Nitromethane	5.3
B3	chloroform	3.9
B4	1,2-dimethoxyethane	4.9
B5	trichloroethylene	5.3
B6	Pyridine	3.8
B7	2-hexanone	1.0
B8	Tetralin	3.6

Nexis is a trademark of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.



Shimadzu Corporation  
[www.shimadzu.com/an/](http://www.shimadzu.com/an/)

SHIMADZU (Asia Pacific) Pte. Ltd,  
[www.shimadzu.com.sg](http://www.shimadzu.com.sg)

For Research Use Only. Not for use in diagnostic procedure.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country. The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. See <http://www.shimadzu.com/about/trademarks/index.html> for details.

Third party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

04-AD-0256-EN

First Edition: Feb. 2022