

CoreFocus Report

No.366


LC Reversed-phase

Shim-pack Series

Shim-pack[™] XR-ODS

High speed analysis of DNPH-Aldehydes

Keywords: Derivatization, DNPH-Aldehydes

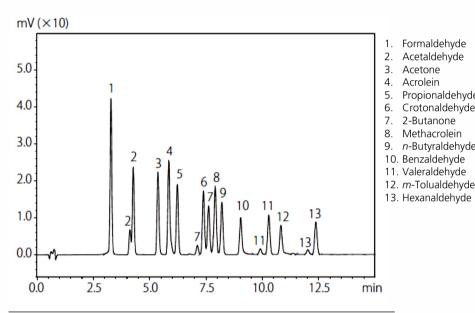

System Prominence™-i (LC-2030C 3D Plus)

Column Shim-pack XR-ODS (75 mm L x 4.6 mm I.D., 2.2 µm)

P/N: 228-41607-93 A) Water/THF = 8/2 (v/v)

B) Acetonitrile

Time	Unit	Parameter	Value
0	Pump	B.Conc	20
15.00	Pump	B.Conc	60
15.01	Pump	B.Conc	20
17.00	Controller	Stop	

Flow rate 1.5 mL/min 40 °C Column temp. Detection UV 360 nm Injection vol. 10 μL

Mobile phase

CoreFocus, Shim-pack and Prominence are trademarks of Shimadzu Corporation or its affiliated companies in Japan and/or other countries

Shimadzu Corporation www.shimadzu.com/an/

For Research Use Only, Not for use in diagnostic procedures,

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®".

Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

Formaldehyde Acetaldehyde

Crotonaldehyde 2-Butanone Methacrolein

n-Butyraldehyde

Acetone Acrolein Propionaldehyde