

# High Capacity Moisture Trap

**1/8-inch tube compression fittings—cat.# 21997**

**1/4-inch tube compression fittings—cat.# 20638**



**Maximum Operating Pressure:** 250 psi

**Maximum Operating Temperature:** 100 °C

## Installation Instructions

1. Remove the protective plug from one end of the trap.
2. Locate nut and ferrule sets (2 each, included). Insert tubing through nut and ferrule until tubing rests firmly against fitting shoulder. Tighten finger-tight.
3. Tighten nut  $3/4$ -turn for  $1/8$ -inch tubing and one full turn for  $1/4$ -inch tubing.
4. Repeat steps 1 through 3 on the other end of the trap.
5. Pressurize the system and leak check with Restek's Electronic Leak Detector (cat.# 22839.)  
DO NOT use liquid leak detectors.

## Health Hazards

**Inhalation:** Slightly irritating to the respiratory tract.

**Skin Contact:** Slightly irritating to the skin.

**Eye Contact:** Slightly irritating to the eyes. May cause redness and pain.

## First-Aid Measures

**Inhalation:** Remove to fresh air.

**Skin:** Wash with soap and water. If irritation persists, contact a physician.

**Eyes:** Flush eyes with plenty of water. Seek medical attention.

**Note:** This document has been prepared based on a component or components of this product. Exposure to the component(s) is not expected through routine laboratory use.

**Call Technical Service at 800-356-1688 or 814-353-1300, ext. 4 (or your Restek representative) if you have any questions about this product or any other Restek product.**

**RESTEK**

Chromatography Products  
© 2010 Restek Corporation

ISO 9001:2008  
cert. # FM80907

#204-08 [002]  
Rev. date: 7/10

**Restek U.S.** • 110 Benner Circle • Bellefonte, PA 16823  
phone: 814-353-1300 • 800-356-1688 • fax: 814-353-1309 • [www.restek.com](http://www.restek.com)

**Restek France** • phone: +33 (0)1 60 78 32 10 • e-mail: [restek@restekfrance.fr](mailto:restek@restekfrance.fr)

**Restek GmbH** • phone: +49 (0)6172 2797 0 • e-mail: [info@restekgmbh.de](mailto:info@restekgmbh.de)

**Restek Ireland** • phone: +44 (0)2890 814576 • e-mail: [restekurope@aol.com](mailto:restekurope@aol.com)

**Restek Japan** • phone: +81 (3)6459 0025 • e-mail: [ryosei.kanaguchi@restek.com](mailto:ryosei.kanaguchi@restek.com)

**Thames Restek U.K. LTD** • phone: +44 (0)1494 563377 • e-mail: [sales@thamesrestek.co.uk](mailto:sales@thamesrestek.co.uk)