



Sample Preparation for GC-MS

Dipl.- Ing. Petra Gerhards
Regional Marketing Manager Columns and Consumables, EMEA
Thermo Fisher Scientific, Dreieich/Germany


Sample Preparation

- Objectives of sample preparation
- Types of sample preparation
 - Fundamentals for different sorbents
 - Polymeric Sorbents
 - Silica based Sorbents
- Fundamentals of solid phase extraction (SPE)
- Solid phase extraction products
 - Protocols
 - Real examples


Not for use in diagnostic procedures

Objectives of Sample Preparation


Concentrate
the analyte

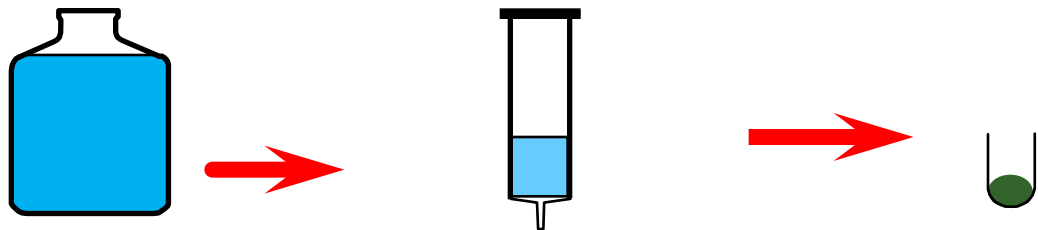

Clean up the
sample matrix

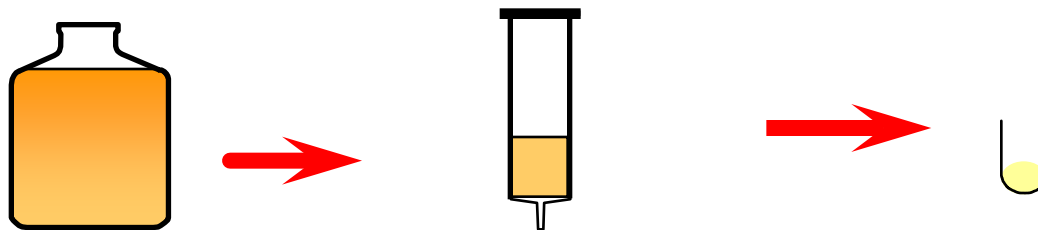

Analytical
compatibility


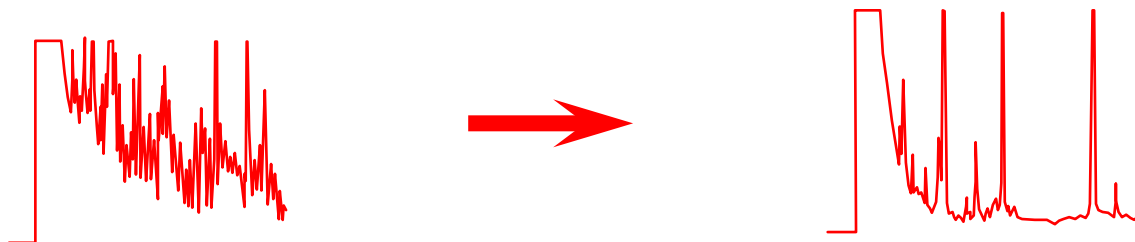


Not for use in diagnostic procedures

Liquid/Liquid Extraction (LLE)

- Does not concentrate the analyte
- Provides a cleaner extract than precipitation
- Analytical compatibility is questionable
 - Note every drug is extracted with the given/chosen solvent for LLE
 - The solvent is always for a large variety of drugs, special drugs like THC require special sample preparation


Analytical
compatibility

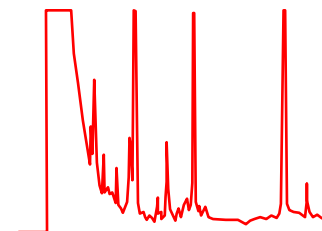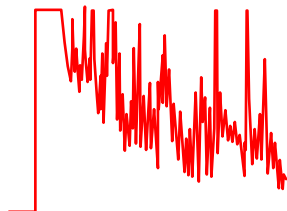

Not for use in diagnostic procedures

Liquid/Liquid Extraction (LLE) with Chlorobutane - Comparison

LLE extracts a lot of hormones and steroids, SPE is MUCH cleaner!

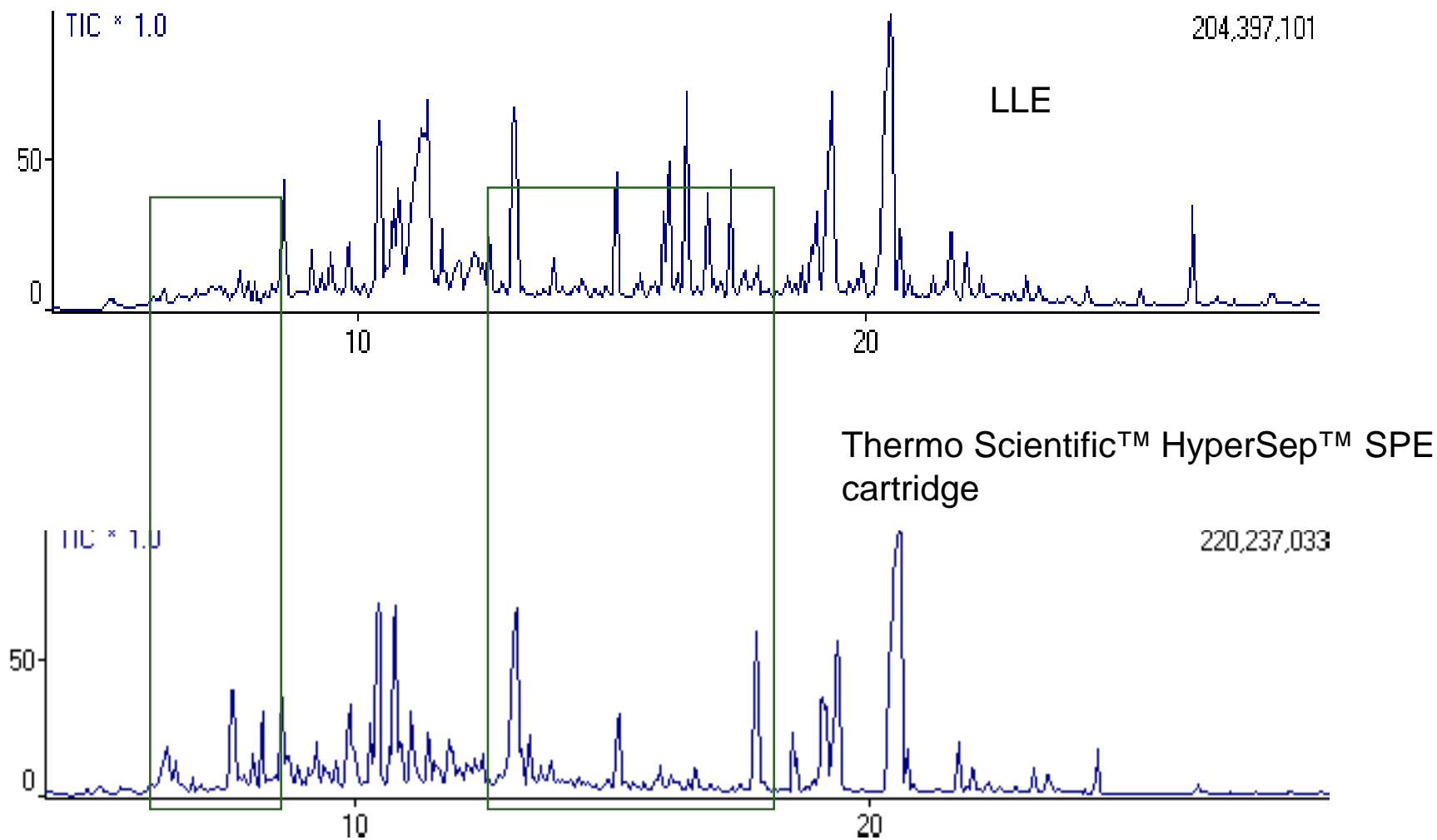

Not for use in diagnostic procedures


Solide Phase Extraction (SPE)

- Capable of concentrating the analytes
 - Up to 2ml of plasma or 3 ml of urine can be used, compared to a 1:1 extraction with LLE
 - Concentration factor after SPE process is 1:60
- Yields very clean extracts
- Extracts are analytically compatible

Not for use in diagnostic procedures

SPE Fundamentals

- Analytes, sample Matrix and SPE sorbent
- Chemistry of extraction protocols
- Troubleshooting and technical comments

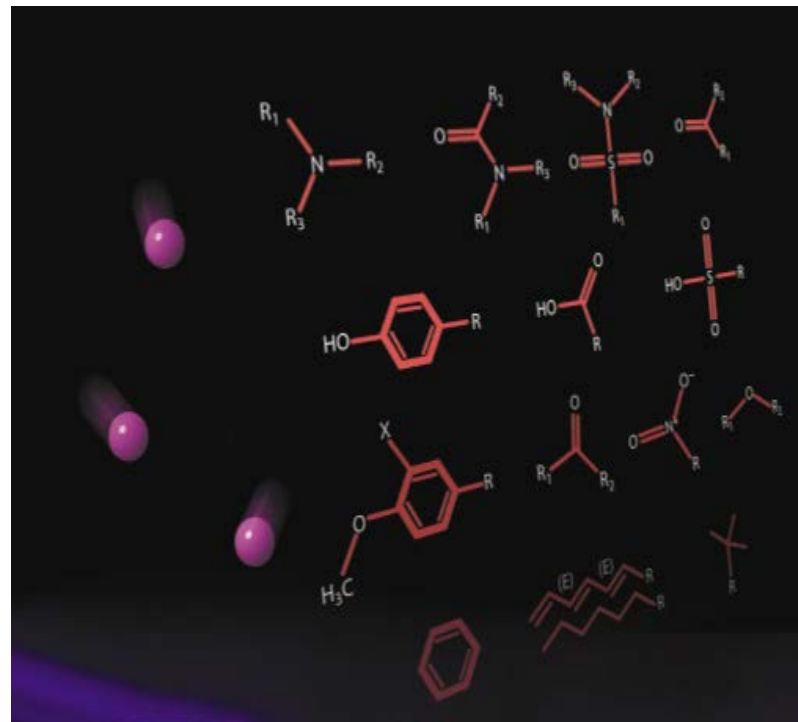

Not for use in diagnostic procedures

Extraction modes

- Nonpolar
- Polar
- Ion exchange
- Covalent*
- Affinity*

Format considerations

- Sample size
- Processing speed
- Device capacity

* Applies to liquid chromatography


Not for use in diagnostic procedures

How Does Chemistry in SPE Work?


$$\text{Log SP} = c + mV_x + rR_2 + s\pi_2^H + a\Sigma\alpha_2^H + b\Sigma\beta_2^H$$




H-Bond Donor

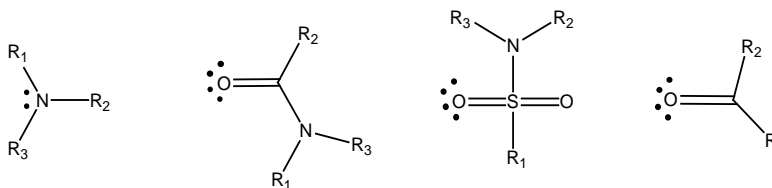


H-Bond Acceptor

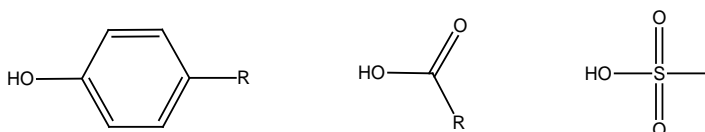

Dipole

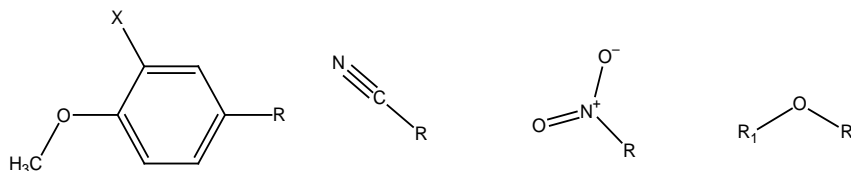


Hydrophobic

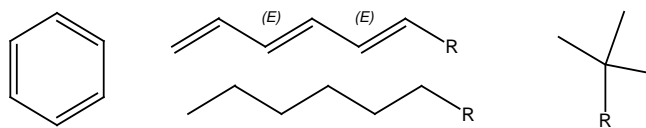

Not for use in diagnostic procedures

Analyte Fundamentals


- Polarity
- Metabolism
- Concentration
- Stability

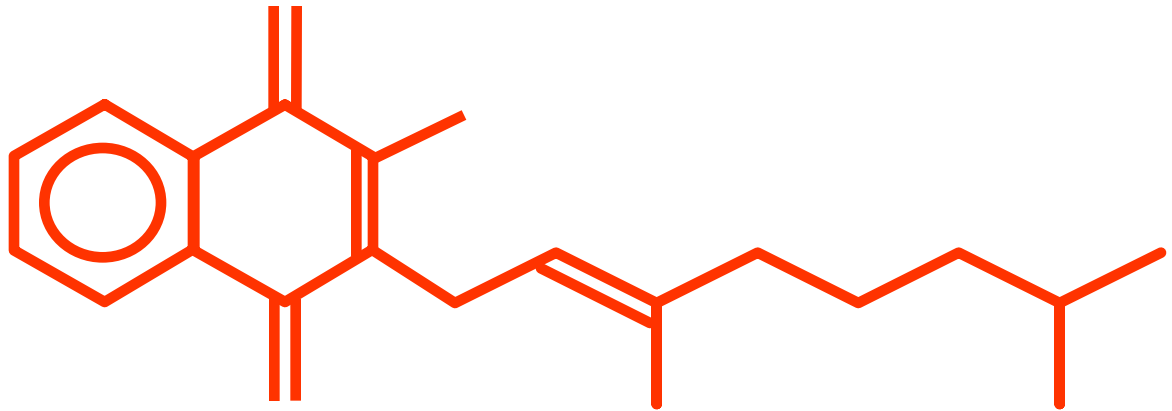

Not for use in diagnostic procedures

Concentration

- What is the maximum concentration expected?
- What is the capacity of sorbent?
- Are there any matrix interferences?

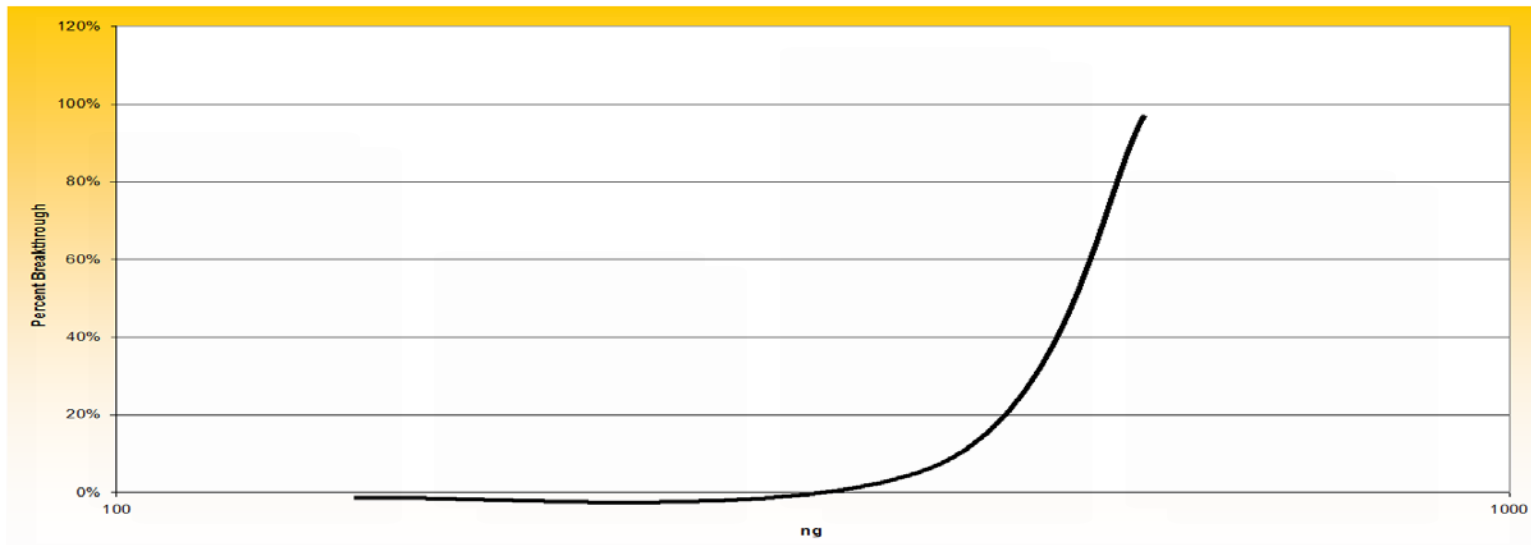

Not for use in diagnostic procedures

Sorbent Capacity

- Capacity is dependent on sorbent type, bed volume and chemistry
- General rule is 4 – 10% of the bed volume on non-polar chemistries
- Certificates of analysis generally indicate sorbent capacity

Not for use in diagnostic procedures

Thermo Scientific™ HyperSep™ and SOLA™ SPE Sorbents

- Silica based → HyperSep cartridges
- Polymeric based → SOLA cartridges

$$\text{Log SP} = c + mV_x + rR_2 + s\pi_2^H + a\Sigma\alpha_2^H + b\Sigma\beta_2^H$$

Not for use in diagnostic procedures

Silica Based Sorbents


- Non-polar (Hydrophobic) phases
 - C18(ec), C18, C8, C4 etc.
- Polar and ion-exchangers
 - CN, Phenyl, SAX, NH₂, SCX, SI
- Mixed phases
 - C8 / SCX



Not for use in diagnostic procedures

Troubleshooting

- Sample pre-treatment → Right pH?
- Conditioning → Is the flow slow enough to activate the sorbent?
- Sample application → pH correct, flow correct?
- Washing → Break through?
- Elution → Solvents correct, enough elution steps?

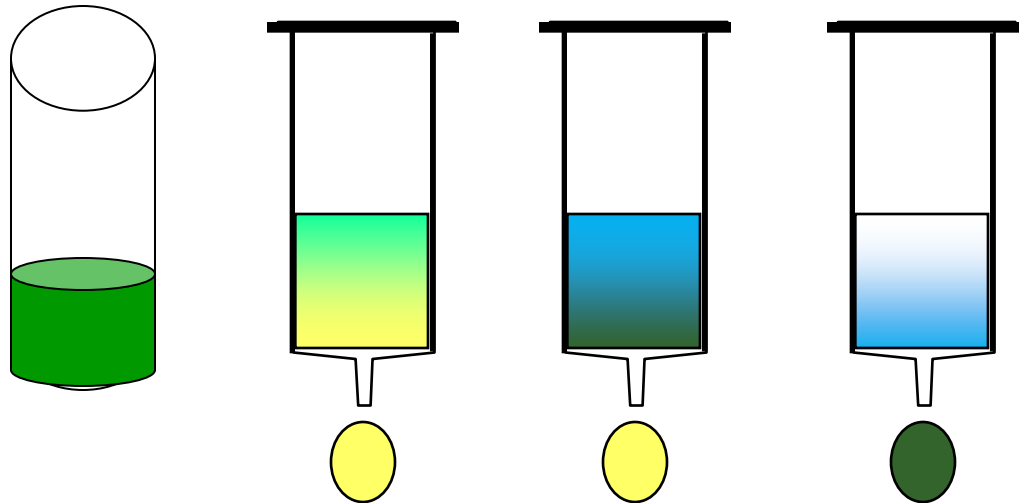

Not for use in diagnostic procedures

HyperSep Based On Cation Exchanger Mechanism


If the substance contains NH or NH₂ or OH groups

- Recommend HyperSep Verify, SCX or WCX will work for these functional groups
- If the substance is a little bit more polar (containing OH groups), recommend SCX
- If the substance is a little bit more non polar, recommend WCX


For ion exchange mechanisms the target is to charge the functional groups.

Not for use in diagnostic procedures


Basic Drugs, Where Charges Can Be Applied

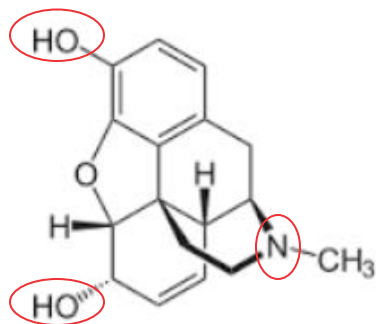

Morphine → SCX or Verify

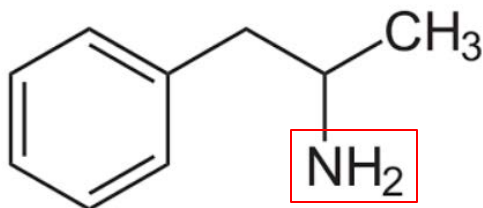

Amphetamine → WCX or Verify

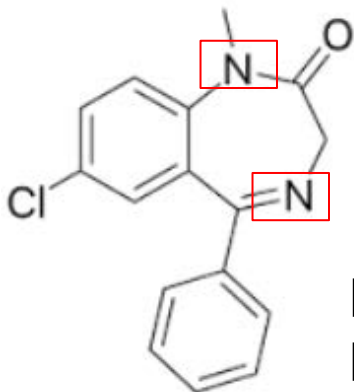


Diazepam → WCX or Verify

For a broad extraction of Drugs of Abuse → **Verify**
For a specific extraction of groups → **SCX or WCX**

Not for use in diagnostic procedures

Zwitterions

If a compound contains NH or NH₂ plus CO, OH or COOH groups recommend HyperSep WCX if the NH or NH₂ group is not steric hindered.

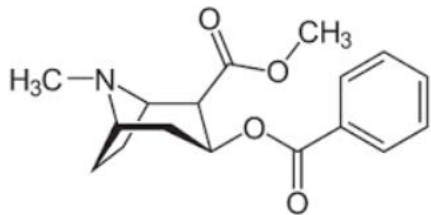


Steric hindered N in Ecgonine
Alakloide from Cocain and
Benzoylecgonine, main metabolite

If it is steric hindered, use HyperSep C18 or HyperSep C18 acid wash, C8 can also be an option.

A method development set would be a good choice, containing HyperSep C18, C8 and Verify.


Not for use in diagnostic procedures

Acidic Drugs

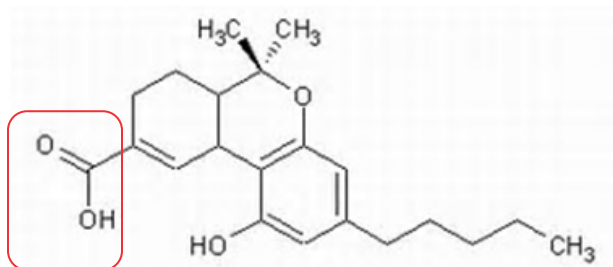

THC carboxylic acid
More hydrophobic interaction

Mostly hydrophobic → C18
For GC sample has to be evaporated
and derivetized and reconstitution
is done on Hexan: Ethylacetate 75:25

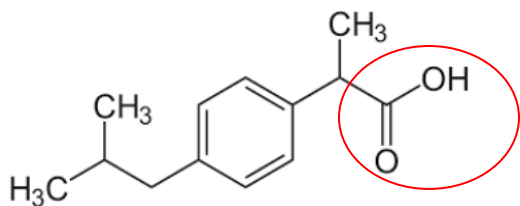

Ibuprofen, acidic drug → **C18 or Verify AX**

Not for use in diagnostic procedures

Cation Exchanger SOLA 10mg/1ml For Plasma Generic Protocol

Low volumes only!


200 µl plasma dilt.

1:1 with dist. water,
add 20 µl acetic acid, pH 4

Interference-
elution

Elution


Condition

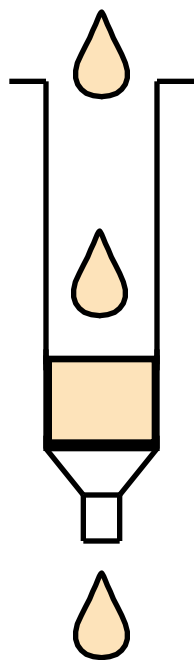

400 µl Methanol
let it react for 2 min.

Equili. with


400 µl 0.1 M acetic acid

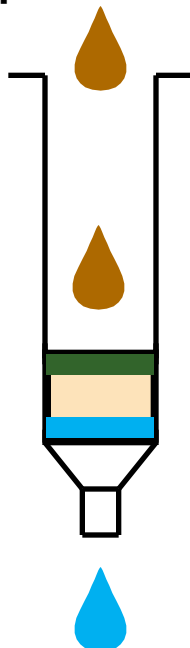

400 µl 0.1 M acetic acid

200 µl methanol

Dry with strong vacuum 2 min.

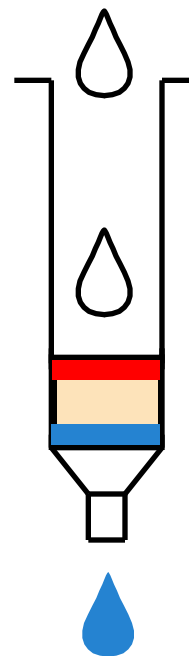

2 x 200 µl


NH₃/methanol

1 : 9

Not for use in diagnostic procedures

SOLA HRP 10mg/1ml For Plasma THX Example

Condition

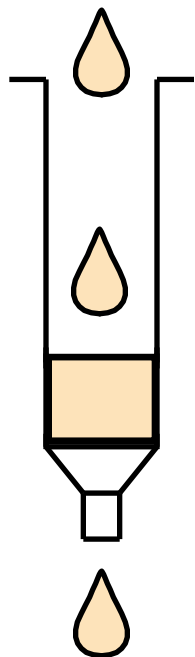


400 μ l Methanol
let it react 2 min.

Followed by 400 μ l neutral
buffer

Ammonia acetate, in some
cases water will be enough

200 μ l plasma dilt.
1:1 with buffer or dist.
Water ,

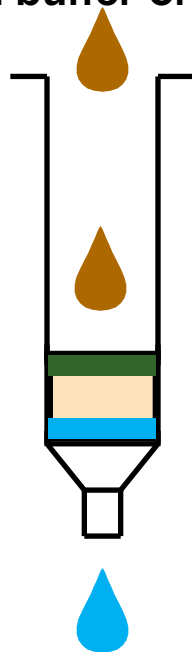


400 μ l buffer + 3 % isopropanol

400 μ l buffer

Dry with strong vacuum 2 min.

Interference-
elution

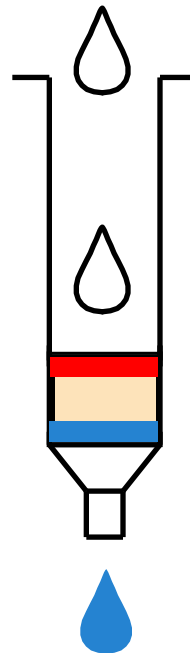

Elution



2 x 200 μ l


Buffer * + ACN

Evaporate, derivetize
and reconstitute in

Hexan 75:Ethylacetate 25


Not for use in diagnostic procedures

THC Sample After Derivatization

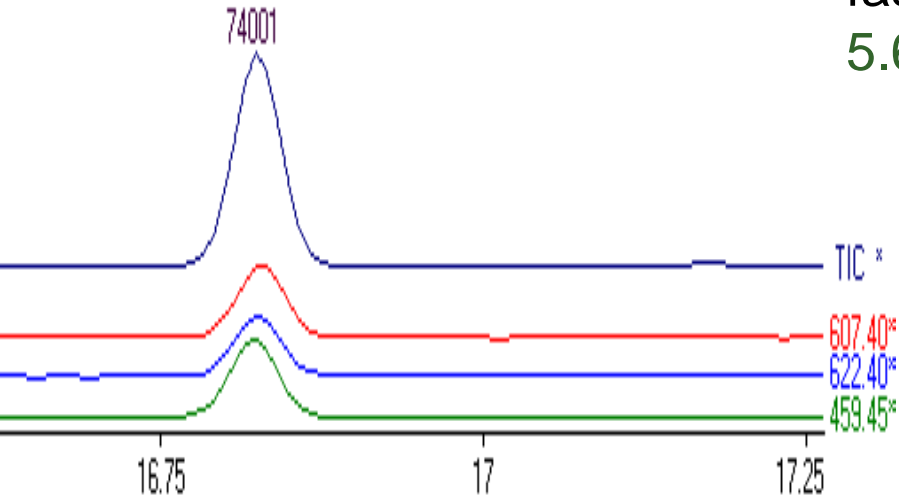

Urine sample
normal GC/MS
16.78 min

Deuterated standard
fast GC/MS
5.68 min

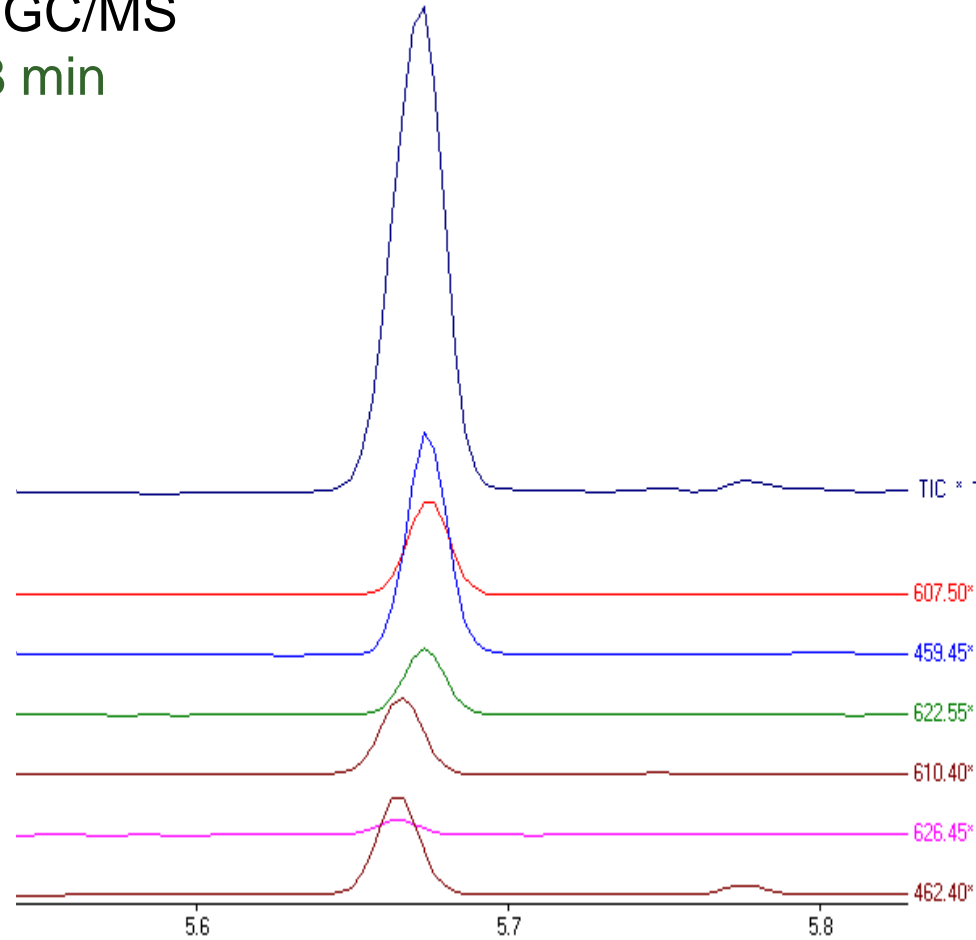

Not for use in diagnostic procedures

HyperSep Verify 200 mg For Urine Generic Protocol

3 ml urine +30 µl

β-Glucuronidase


Incubate at 56 ° C for 60 min

Analyte elution with

2 x 2 ml

NH₃/methanol

Sample application

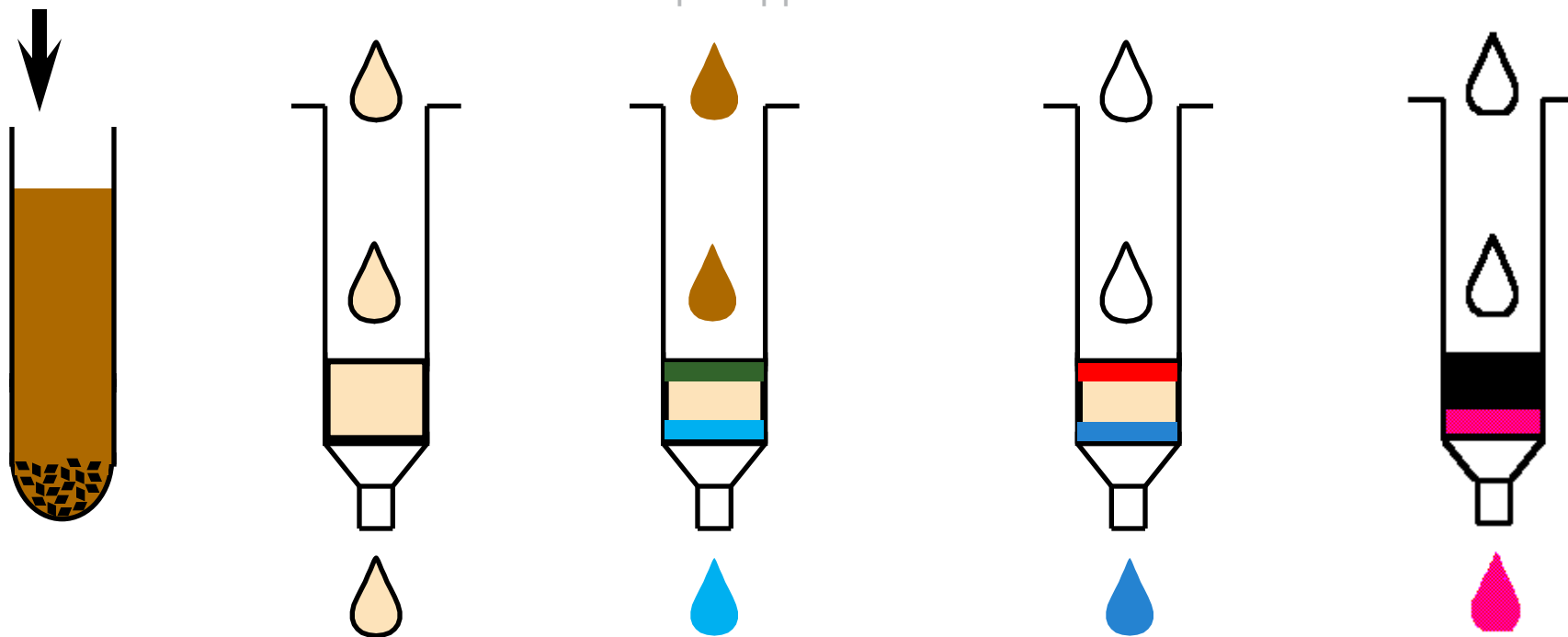

Conditioning with 3ml methanol
pre-equilibration with
diluted phosphoric acid

Interference elution with 3 ml

Diluted phosphoric acid, 0.5 ml 0.1 N acidic acid and


1ml methanol

Not for use in diagnostic procedures

SPICE Urine WITH ME OF Synthetic Cannabinoides

3.8. 11.20 Uhr Crea 45.2 mM/l
3.8. 18.45 Uhr Crea 3.6 mM/l
4.8. 10.00 Uhr Crea 14.7 mM/l

Without SPE you can not determine
SPICE ME, concentration is too small,
within 24 hrs the ME are already excreeded

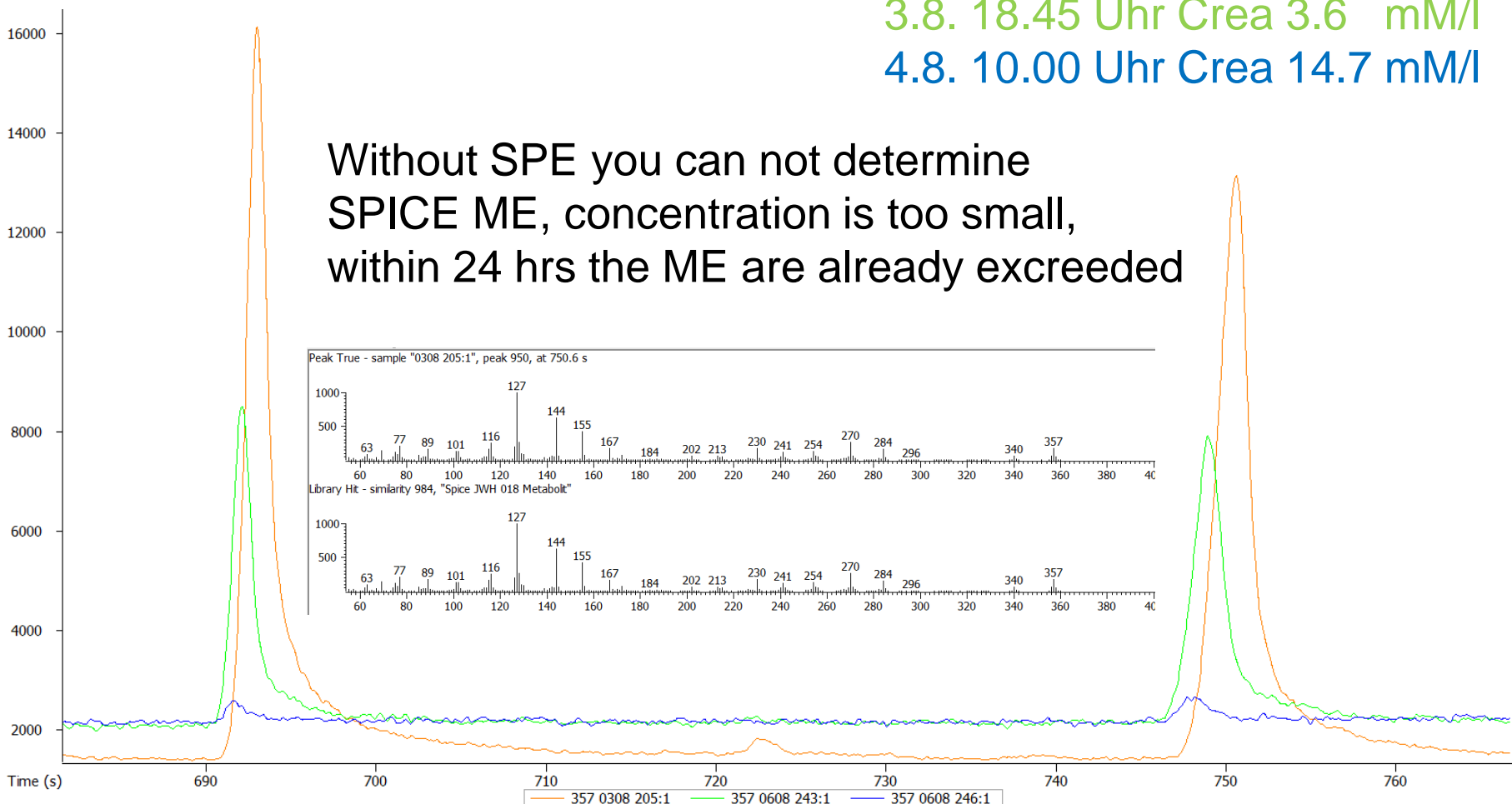

Not for use in diagnostic procedures

Any Questions?

Do you want to receive detailed SPE protocols?
We have protocols ready for drugs of abuse analysis.
Others could be assembled upon request.
Please send an e-mail to
petra.gerhards@thermofisher.com and specify
if required for GC or LC.



**Do you have additional questions
or do you want to talk to an expert from Thermo Fisher
Scientific?**
Please send an E-Mail to analyze.eu@thermofisher.com
and we will get back to you.